

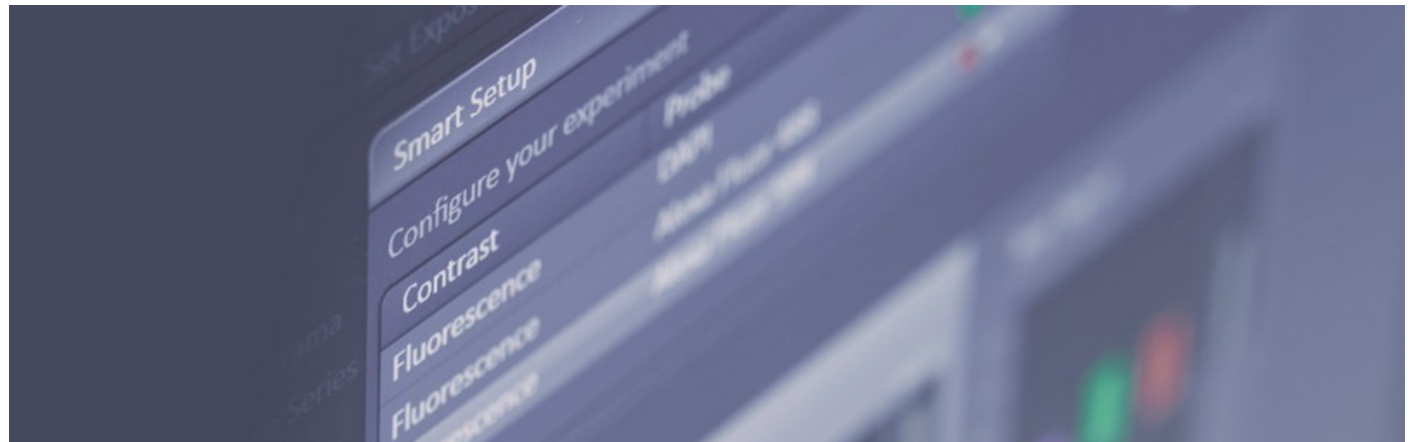


How to Create the Light Source Short Cut Buttons

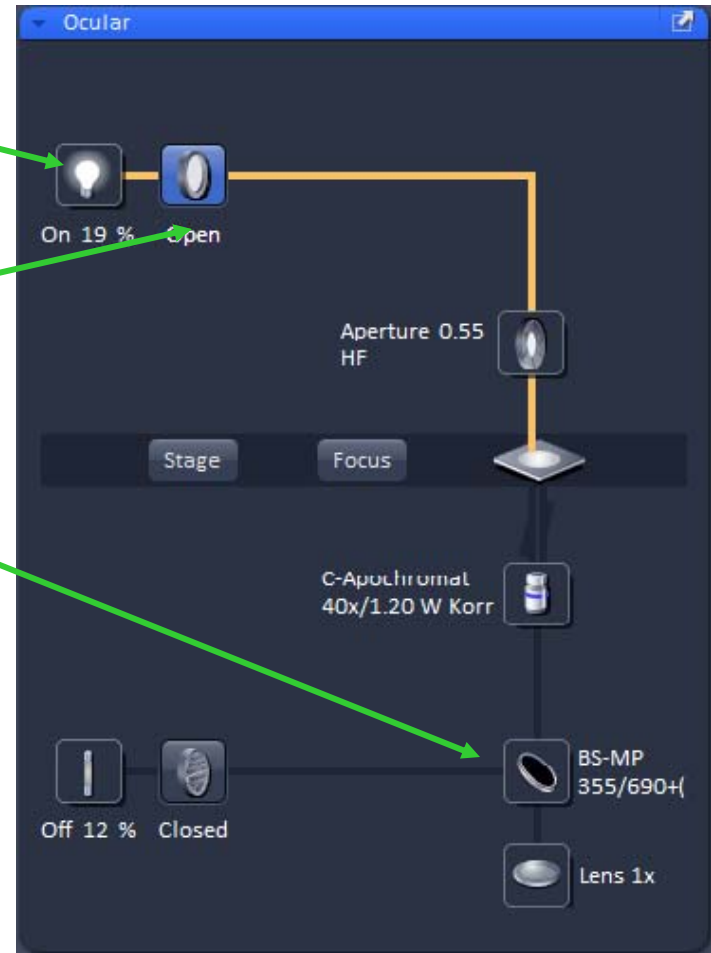


Transmitted Light ON

- Select 'Oculars online'
- Turn on the transmitted light source.
- Open the transmitted light shutter.
- Select BS-MP 355/690
- Save the configuration

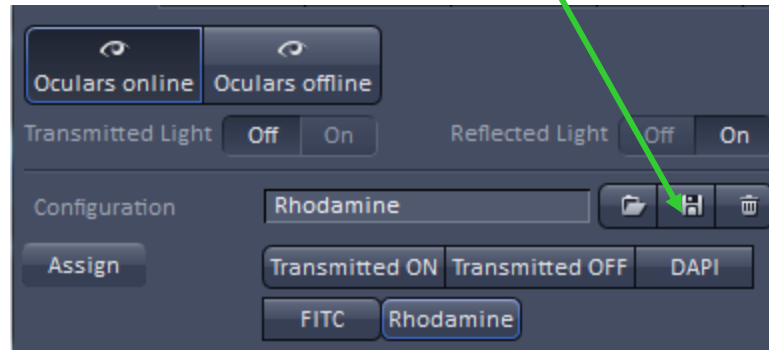



Transmitted Light OFF

- Turn off the transmitted light source.
- Close the transmitted light shutter.
- Save the configuration

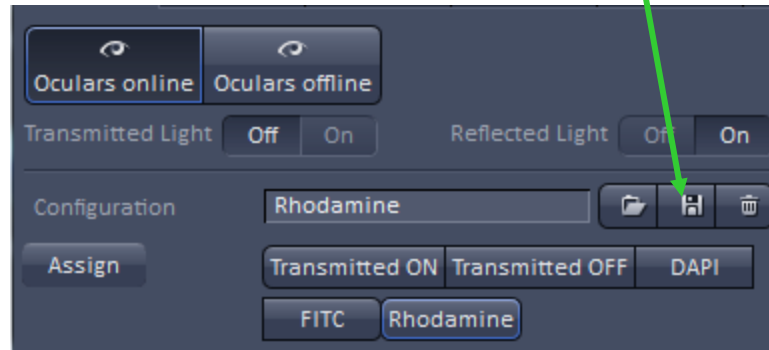



DAPI Light Setting

- Turn on the fluorescent light source, increase the power to 12%.
- Open the fluorescent light shutter.
- Select the DAPI filter.
- Save the configuration.

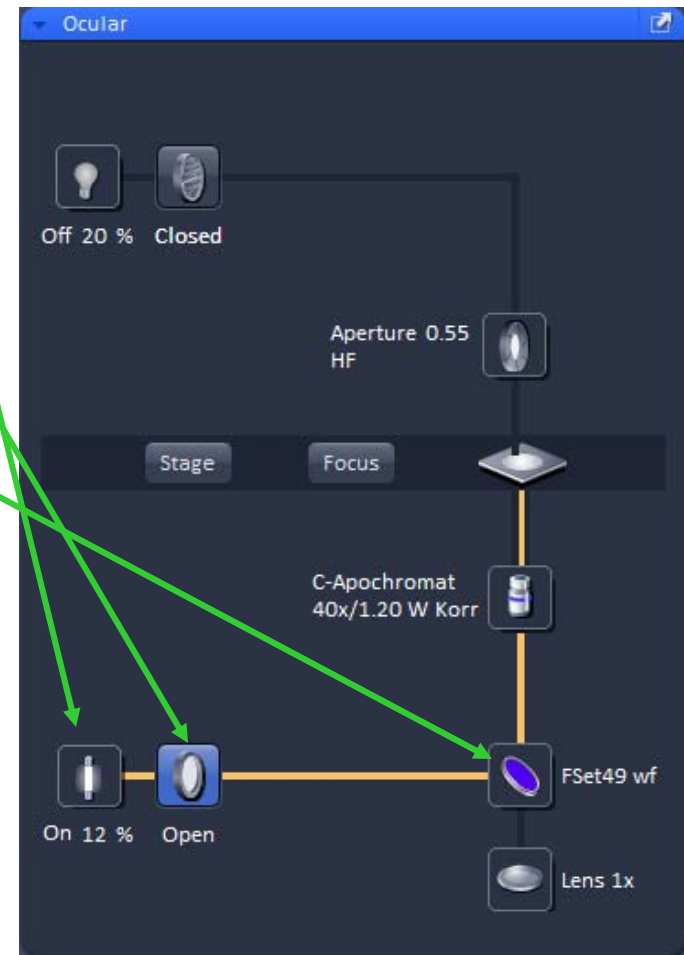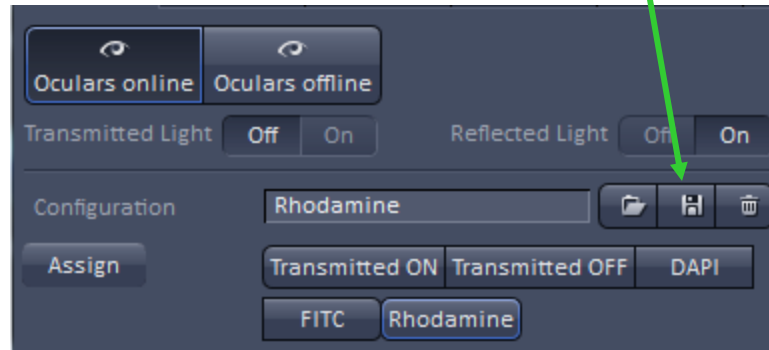



Green Light Setting

- Turn on the fluorescent light source, increase the power to 12%.
- Open the fluorescent light shutter.
- Select the Green filter.
- Save the configuration.

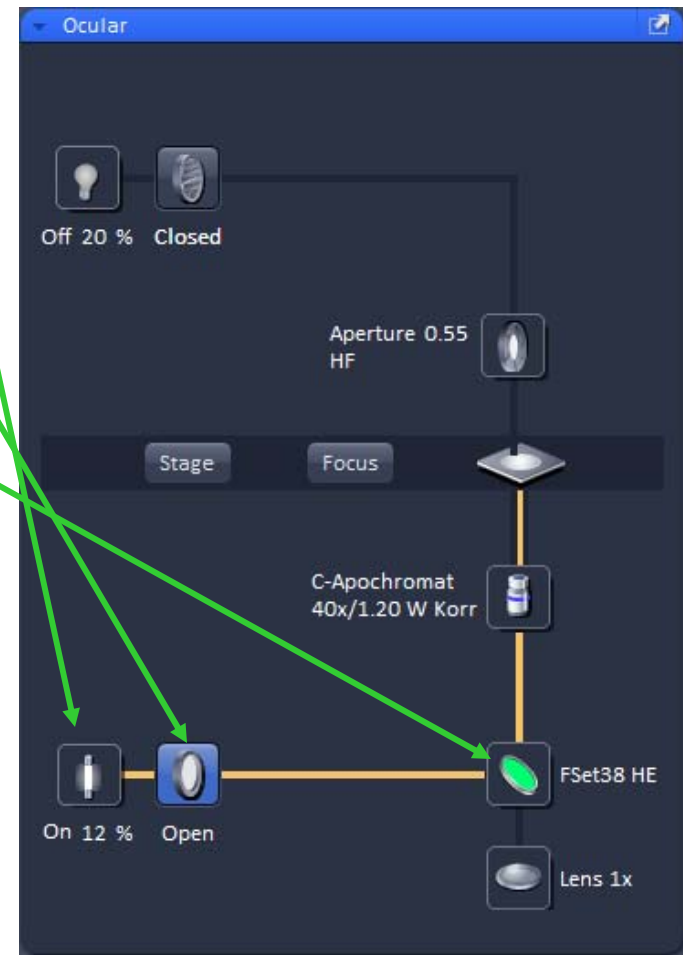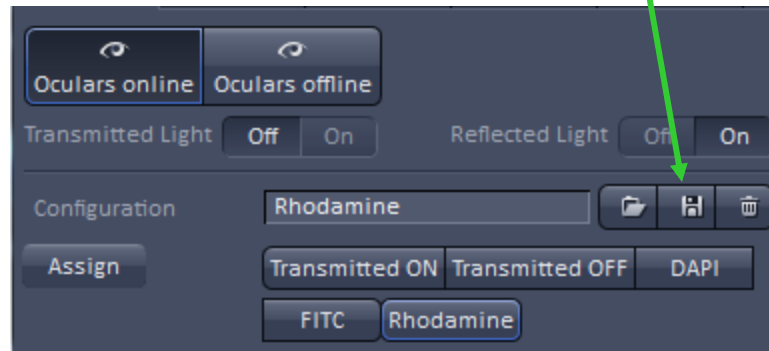

Red Light Setting

- Turn on the fluorescent light source, increase the power to 12%.
- Open the fluorescent light shutter.
- Select the Red filter.
- Save the configuration.

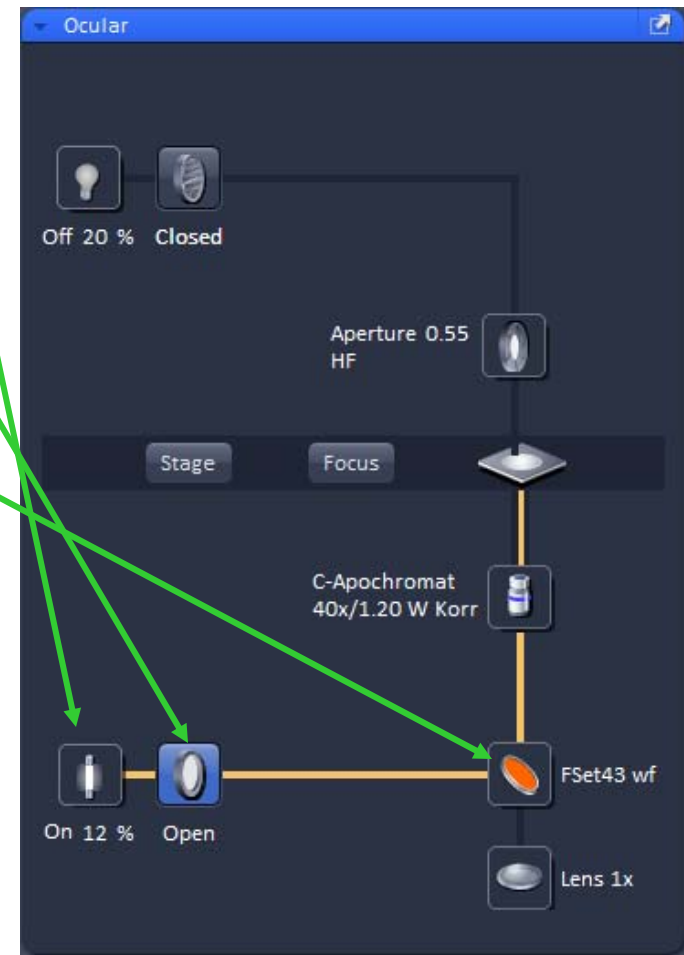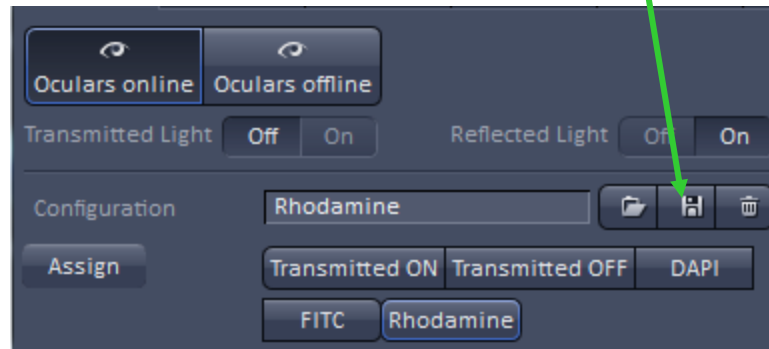

Fluorescent Light OFF/Creating Buttons

- Turn off the fluorescent light source, for each color, by pressing 'OFF'.
- Select 'Assign'
- Select Button #
- Select Configuration
- Select Assign
- Perform for Transmitted Light ON and OFF, DAPI, Green and Red filter configurations.

